Product code: smsG1B-1




Introduction

The ChromoTek Nano-CaptureLigand™ mouse IgG1, Fc-specific VHH, biotinylated is used for the site-directed and specific immobilization of the Fc-fragment of mouse IgG1 in biosensor and ELISA assays. It captures non-biotinylated mouse IgG1 antibodies or Fc-fragments to streptavidin/avidin.

Nano-CaptureLigand mouse IgG1, Fc-specific VHH, biotinylated comprises a monoclonal biotinylated VHH/ Nanobody. The product belongs to the Nano-CaptureLigands™ family.

Properties

Description	Monovalent, recombinant single domain antibody for the immobilization of mouse IgG1: alpaca monoclonal Nanobody, Fc-specific, biotinylated	
Product Type	Capture Nanobody (VHH)	
Applications	Immobilization of mouse IgG1 antibodies on avidin and streptavidin surfaces for Bio-Layer Interferometry (BLI), Surface Plasmon Resonance (SPR) and ELISA	
Target / Specificity	Fc-fragment of mouse IgG1	
Cross-reactivity	No cross-reactivity to mouse IgG2a, Ig2b, IG2c, IgG3; human IgG1, IgG2, IgG3, IgG4, IgA, IgM, IgE; rabbit IgG; rat IgG1, IgG2a, IgG2b, IgG2c; sera from goat, guinea pig, human, macaque (cynomolgus monkey), rabbit, rat, sheep	
Affinity (<i>Kd</i>) of monovalent (1:1) binding mode	0.13 nM Apparent affinity may be higher for full IgGs due to avidity effects (1 antibody captured by 2 Nanobodies).	
Concentration	1 g/L (73 μM)	
Conjugate	Biotin	
Degree of biotinylation	On average 1-2 biotin molecules per Nanobody	
Format	Alpaca single domain antibody, monovalent	
Host	Alpaca-derived, recombinantly produced in bacteria	
Clonality	Monoclonal	
Clone	CTK0103 (VHH0295)	
RRID	AB_2848186	

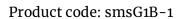



Synonyms	Alpaca single domain antibody, VHH, Nanobody, binding domain of single domain antibody, Nano-antibody
Validation	Application validated for ELISA and BLI (FortéBio Octet® systems) Determination of cross-reactivity, subclass specificity, sequence, affinity, and melting temperature
Purity	Recombinantly expressed and purified via His-tag
Form	Buffered aqueous solution
Storage buffer	25 mM TAPS pH 8.5, 500 mM NaCl, 5 mM EDTA, Preservative: 0.09 % sodium azide
Storage conditions	Upon receipt store at +4°C/+40°F. <i>Optional</i> : Aliquot upon arrival and store at -20°C/-4°F
Stability	Stable for 1 year at +4°C/+40°F
Shipment	Shipped at ambient temperature

Product sizes

Product	Product code	Size
Nana Cantural impadim popusa IgC1. Es apacific VIIII, histiguiated	smsG1B-1-10	10 μL
Nano-CaptureLigand™ mouse IgG1, Fc-specific VHH, biotinylated	smsG1B-1-100	100 μL

Product code: smsG1B-1



Suggested buffer compositions

Recommended buffers for BLI

Buffer	Composition	
1x Kinetics buffer	PBS, 0.01 % (m/v) BSA, 0.002% (v/v) Tween-20	
Regeneration buffer	0.01 M glycine, pH 2	

Recommended buffers for ELISA

Buffer	Composition	
Blocking buffer	PBS, 3 % BSA	
Dilution buffer	PBS, 0.05 % Tween-20, 0.5 % BSA	
Wash buffer for avidin coating	PBS	
Wash buffer	PBS, 0.05 % Tween-20	

Product code: smsG1B-1



Bio-Layer Interferometry (BLI) protocol

General notes

- The following protocol was designed for the use with a FortéBio Octet Red96e system. Other BLI systems or your specific research question may require optimization of particular parameters.
- Use the recommended materials or their equivalents.
- Set up all samples in a black 96-well microplate (e.g. Greiner Microplate 96 well, PP, flat-bottom, black, #655209) at room temperature. Use 200 μL per well.
- Nano-CaptureLigands are highly compatible with avidin or streptavidin sensors (e.g. FortéBio Streptavidin (SA) Biosensors, #18-5019) and FortéBio Octet[®] and and BLltz[®] systems.
- Run all experiments at +30°C, a shaking speed of 1000 rpm and a recording rate of 5 Hz.
- Dilute all samples in 1x Kinetics buffer. Optional: Use 10x Kinetics buffer.
- NanoCaptureLigands can be regenerated at least 10 times with Regeneration buffer with minimal loss of binding efficiency.
- Nano-CaptureLigands carry a His-tag; thus, avoid the use of anti-His primary antibodies.
- Briefly centrifuge the Nano-CaptureLigand solution before use.

Protocol

1. Baseline 1:

• Incubate the biosensors for 60 s in 1x Kinetics buffer.

2. Loading

- Dilute the Nano-CaptureLigand to a concentration of 1 µg/mL in 200 µL 1x Kinetics buffer.
- Load the diluted Nano-CaptureLigand onto the biosensors for 60-120 s until a loading response of 1 nm is reached.

Optional: Use the threshold limit function in the FortéBio Data Acquisition software.

3. Quenching (optional)

• Incubate the biosensors for 60 s with biocytin (10 µg/mL in 1x Kinetics buffer).

4. Baseline 2

• Incubate the biosensors for 120 s in 1x Kinetics buffer.

5. Activation

• Activate the biosensors for 120-180 s with the antibody (20 nM in 1x Kinetics buffer).

Product code: smsG1B-1



6. Baseline 3

• Incubate the biosensors for 120 s in 1x Kinetics buffer.

7. Association

• Bind different antigen concentrations in 1x Kinetics buffer for 120-600 s. *Note:* As a start, use 0.1-250 µg/mL or 1/10-10x *Kd* of antigen.

8. Dissociation

• Incubate the biosensors for 60-800 s in 1x Kinetics buffer.


Note: Use the wells from step 6, Baseline 3.

Note: Duration of dissociation step depends on the affinity of the analyzed interaction.

9. Regeneration (optional)

- Regenerate the biosensors for 5 s with Regeneration buffer.
- Incubate in 1x Kinetics buffer for neutralization.
- Repeat regeneration 2 times.


Application examples

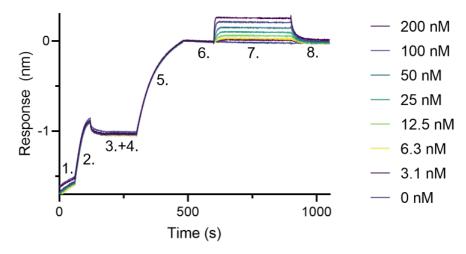

BLI binding kinetics of a mouse IgG1 anti-DYKDDDDK-tag antibody to a DYKDDDDK-tagged protein. A monoclonal mouse IgG1 anti-DYKDDDDK-tag antibody was immobilized using Nano-CaptureLigand mouse IgG1, Fc-specific VHH, biotinylated on FortéBio Streptavidin (SA) Biosensors and assayed with different concentrations of a DYKDDDDK-tagged protein.

Product code: smsG1B-1



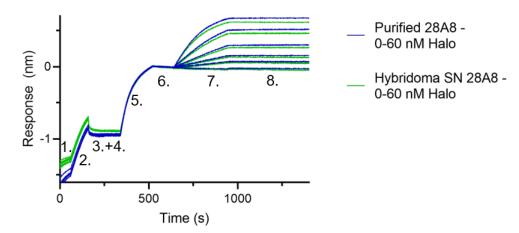

BLI binding kinetics of a purified and a hybridoma supernatant mouse IgG1 anti-Halo antibody to Halo-tag. A purified, monoclonal mouse IgG1 anti-Halo antibody (ChromoTek Halo antibody [28A8]) and a hybridoma supernatant, monoclonal mouse IgG1 anti-Halo antibody were immobilized using Nano-CaptureLigand mouse IgG1, Fc-specific VHH, biotinylated on FortéBio Streptavidin (SA) Biosensors and assayed with different concentrations of Halo-tag protein.

Product code: smsG1B-1



Sandwich ELISA protocol

General notes

- The following protocol was designed for a standard sandwich ELISA. Other types of ELISA or your specific research question may require optimization of particular parameters.
- Use the recommended materials or their equivalents.
- In this protocol, MaxiSorp plates (e.g. Thermo Scientific™ White and Black 384-Well Immuno Plates, #460518) are used that must be coated with avidin or streptavidin first. Alternatively, pre-coated avidin/streptavidin plates can be used.
- Nano-CaptureLigands carry a His-tag; thus, avoid the use of anti-His primary antibodies.
- Recommended volumes for 96-well and 384-well microplates:

Protocol steps	96-well microplate	384 -well microplates
Coating, antigen binding, antibody binding	100 μL	20 μL
Washing, blocking	300 μL	90 μL

• Briefly centrifuge the Nano-CaptureLigand solution before use.

Protocol

1. Avidin coating (optional)

- Coat each well of a MaxiSorp plate with 10 µg/mL avidin in PBS at +4°C overnight.
- Wash each well twice with PBS.

2. Blocking

- Block each well with Blocking buffer for 1-2 h at room temperature.
- Wash each well 3 times with Wash buffer.

3. Nano-CaptureLigand coating

- Add 50 nM Nano-CaptureLigand (diluted in Dilution buffer) to each well.
- Incubate for 1 h at room temperature.
- Wash each well 5 times with Wash buffer.

4. Immobilization of capture antibody

- Add the capture antibody (diluted in Dilution buffer) to each well.
- Incubate for 1 h at room temperature.
- Wash each well 5 times with Wash buffer.

Note: Test different concentrations of the capture antibody in an initial experiment.

Product code: smsG1B-1



5. Antigen binding

- Add the antigen to each well.
- Incubate for 1 h at room temperature.
- Wash each well 5 times with Wash buffer.

Note: Test different concentrations of the antigen.

6. Binding of primary antibody

- Add the primary antibody to each well and incubate.
- Wash each well 5 times with Wash buffer.

Note: Dilute and incubate the primary antibody as indicated in the manufacturer's manual.

7. Binding of secondary / detection antibody

- Add the secondary / detection antibody to each well and incubate.
- Wash each well 5 times with Wash buffer.

Note: Dilute and incubate the secondary / detection antibody as indicated in the manufacturer's manual.


8. Detection

- Add the appropriate ELISA substrate solution to each well and incubate as indicated in the manufacturer's manual.
- Analyze with a microplate reader.

Product code: smsG1B-1



Application examples

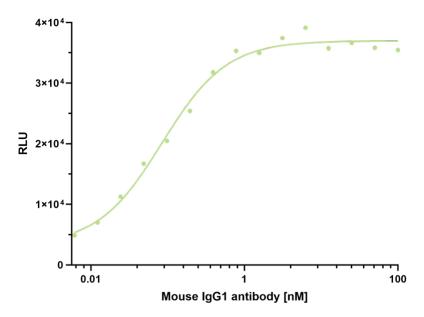

ELISA capture of a mouse IgG1 antibody using Nano-CaptureLigand mouse IgG1, Fc-specific VHH, biotinylated. 50 nM Nano-CaptureLigand mouse IgG1, Fc-specific VHH, biotinylated was used for coating on an avidin-coated MaxiSorp plate. Mouse IgG1 antibody was titrated in a 1:2 dilution series and detected with an alkaline phosphatase-conjugated detection antibody.

Product code: smsG1B-1



Product overview and related products

Product	Product code
Nano-CaptureLigand™ human IgG/rabbit IgG, Fc-specific VHH, biotinylated	shurbGB-1-10; -100
Nano-CaptureLigand™ human Ig, lambda-LC-specific VHH, biotinylated	shuLB-1-10; -100
Nano-CaptureLigand™ human IgE, VHH, biotinylated	shuEB-1-10; -100
Nano-CaptureLigand™ mouse IgG1, Fc-specific VHH, biotinylated	smsG1B-1-10; -100
Nano-CaptureLigand™ mouse IgG2a, Fc-specific VHH, biotinylated	smsG2aB-1-10; -100
Nano-CaptureLigand™ mouse IgG2b, Fc-specific VHH, biotinylated	smsG2bB-1-10; -100
Nano-CaptureLigand™ mouse IgE, VHH, biotinylated	smsEB-1-10; -100
GFP VHH, biotinylated recombinant binding protein	gtb-250

For product details, information, and ordering visit www.chromotek.com.

Product code: smsG1B-1



Contact

support@chromotek.com

ChromoTek GmbH Am Klopferspitz 19 82152 Planegg-Martinsried Germany phone: +49 89 124 148 80 fax: +49 89 124 148 811 ChromoTek Inc. 62-64 Enter Lane Islandia, NY 11749 USA phone: 631 501 1058 fax: 631 501 1060

Disclaimer

Only for research applications, not for diagnostic or therapeutic use!

ChromoTek and GFP-Trap®, RFP-Trap®, Myc-Trap®, Spot-Trap®, Spot-Trap®, Spot-Label®, Spot-Cap®, Nano-Secondary®, F2H® Kit, and Chromobody® are registered trademarks and Nano-CaptureLigand™ is a trademark of ChromoTek GmbH. Nanobody is a registered trademark of Ablynx, a Sanofi company. Octet® is a registered trademark of ForéBio, a Sartorius brand. Dynabeads™ is a trademark of Life Technologies AS, a part of Thermo Fisher Scientific Inc. Alexa Fluor® is a registered trademark of Life Technologies Corporation, a part of Thermo Fisher Scientific Inc. Other suppliers' products may be trademarks or registered trademarks of the corresponding supplier each. Statements on other suppliers' products are given according to our best knowledge.